


Notes: _____

DIGITAL SCALES

MODEL DKS-3002 OPERATING INSTRUCTIONS

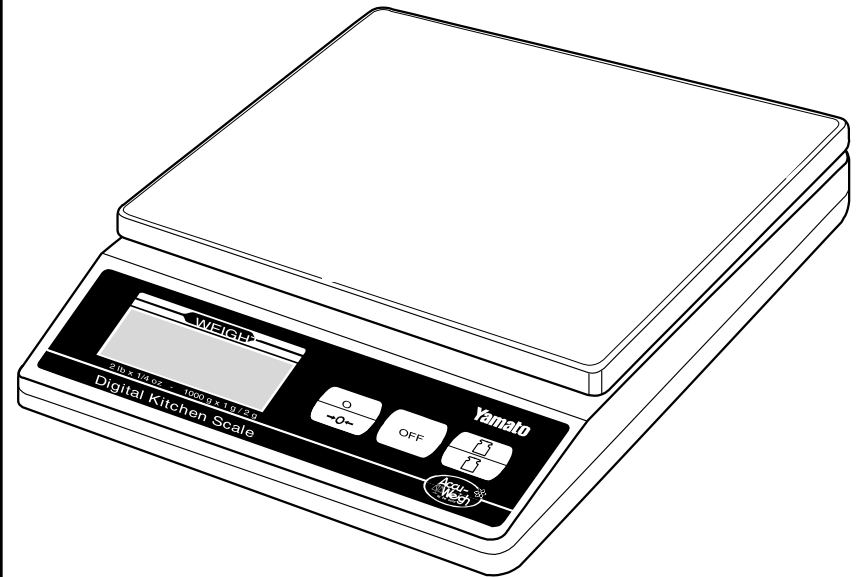

by **Yamato**

Yamato
www.YamatoCorp.com


YAMATO CORPORATION
1775 S. Murray Blvd.
Colorado Springs, CO 80916 USA
Tel (719) 591-1500 Fax (719) 591-1045

YAMATO TECH CORPORATION
#112-19425 Langley By-Pass
Surrey, B.C. V3S 6K1 Canada
Tel (604) 533-2338 Fax (604) 533-0827

INSTRUCTIONS FOR USE


TO FIT BATTERY

Remove the cover on the underside of the scale and push the connector firmly onto the battery terminals. Refit the cover. (Use 9 V Alkaline battery for suitable power output.)

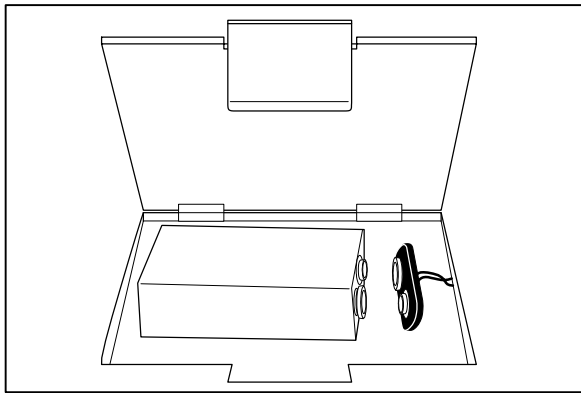

TO SWITCH ON

Place the scale on a flat, stable surface. Press the ON/Zero key once (quite firmly) in the center. The display will show "8888" then "0:0" (or "0".)

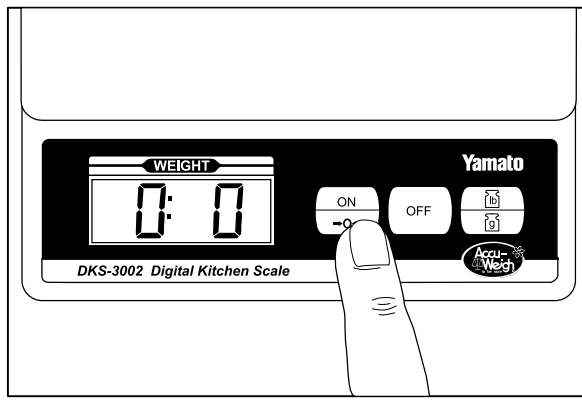

CLEANING AND CARE

- If the scale is not being used for a long period of time, remove the battery. Remove dead batteries immediately.
- Clean scale with damp cloth, do not immerse the scale in water or use chemical/abrasive cleaning agents.
- All plastic parts should be cleaned immediately after contact with fats, spices, vinegar or strongly flavored/colored foods. Avoid contact with acids such as citrus fruit juices.
- Avoid placing your scale directly next to a microwave oven if both will be used at the same time. (Some microwave ovens emit interference which, at close range, can cause the display of the scale to fluctuate.)

TECHNICAL SPECIFICATION

Power Supply:

One 9V Alkaline battery

Accuracy:

To within +/- one increment or 1% of applied load

MINOR TROUBLE SHOOTING

1. Is the scale level?

Make sure that all the feet are resting firmly on the supporting surface.

2. Is the scale turned on?

If the scale does not work at all, depress the ON/Zero key, Repeat once or twice. Scale will automatically turn off if not used for two minutes (thirty seconds with a zero reading) to preserve battery life.

3. "Err" indication?

When the weight of the scale platform exceeds 105% of the capacity of the scale, "Err" will appear on the digital display, Reduce load.

4. Is movement of the scale restricted in any way?



Make sure that no foreign objects are wedged between the platform and the scale body. Verify that platform is not rubbing against surrounding objects.

5. "Lo" indication?

Remove battery cover and install new 9 Volt Alkaline battery. If any of these actions fail to correct the problem, the scale needs to be serviced by a qualified technician.

TO CONVERT WEIGHT UNITS

Press the LB/G key to convert between metric and imperial units. This can be done at any time during weighing.

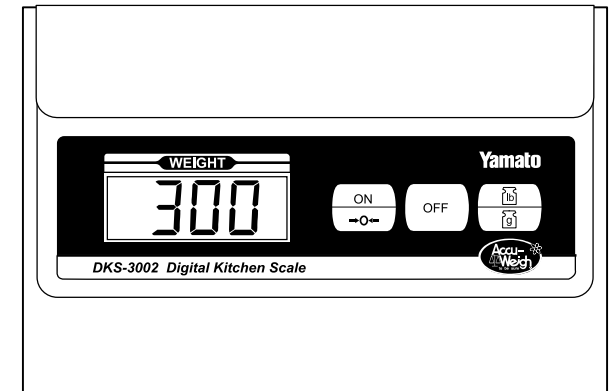
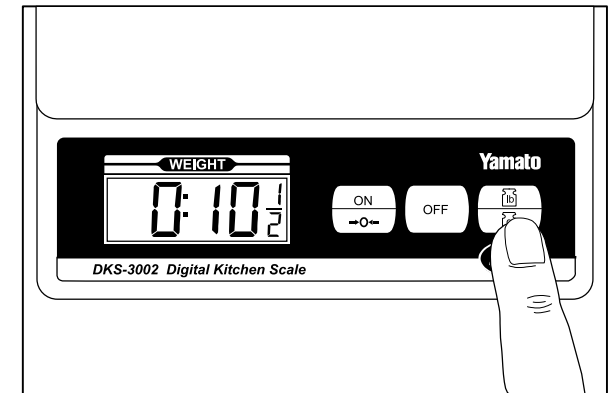


USING A BOWL


This scale may be used with any bowl or container which weighs less than 1 kg or 2 lb 3 oz.

Place the bowl on the platform then press the ON/Zero key.

The display will show "tArE" for 2 seconds, then it will reset to "0:0" (or "0"), ready to weigh.

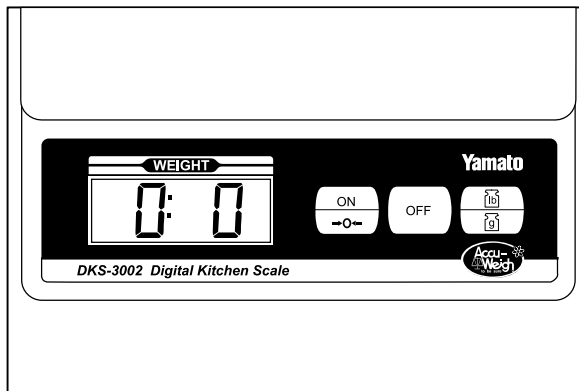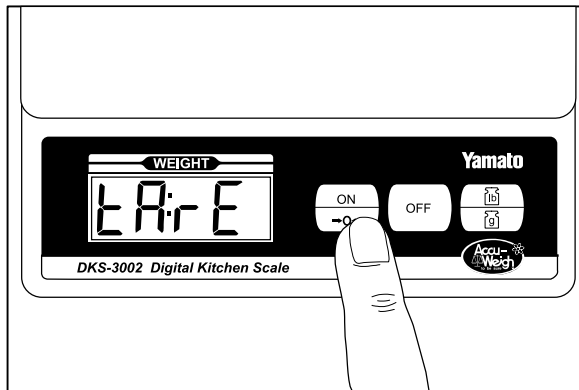**MINUS READINGS**

"----" indicates a minus reading, i.e. less than zero. For example, if the scale is set to zero with a bowl on the platform then the bowl is removed, the display will show "----". To reset to zero, either replace the same bowl or press the ON/Zero key.

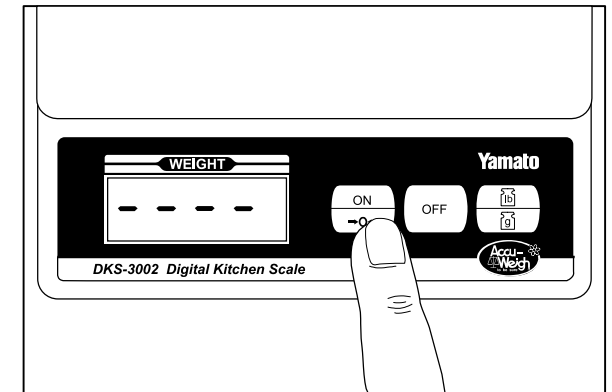**TO SWITCH OFF**

Press the OFF button once to switch off. Switching your scale off immediately after use will increase battery life. An automatic switch-off will operate after 30 seconds if the display has been showing zero continuously for that time, or after 2 minutes if a weight reading has been displayed continuously for that time.

